



# GeneJET™ RNA Purification Kit


Fermentas launched GeneJET™ RNA Purification Kit - a simple and efficient system for purification of total RNA from mammalian cultured cells, tissue, human blood cells, bacteria and yeast.

## Features


- **Universal** – suited for both cell and tissue samples from a wide range of sources
- **Efficient** – high yields of total RNA
- **Fast** – 15 minutes after lysis step
- **Pure** – an  $A_{260/280}$  ratio > 1.9

Product No.	Product Description	Size
K0731	<b>GeneJET™ RNA Purification Kit</b>	50 preps
K0732	<b>GeneJET™ RNA Purification Kit</b>	250 preps


## Overview of Purification Procedure

- 


**Collect sample**  
**Prepare lysate**
- ↓ **Add ethanol**




**Transfer** solution to the GeneJET™ RNA Purification Column

**Centrifuge** 1 min  
**Discard** the flow-through  
**Repeat** step 2
- 

**Transfer** column into a new tube  
**Add** Wash Buffer 1

**Centrifuge** 1 min  
**Discard** the flow-through
- 

**Add** Wash Buffer 2

**Centrifuge** 2 min  
**Discard** the flow-through  
**Repeat** step 4
- 

**Transfer** column into a new tube  
**Add** 50 - 100 µl Water, nuclease-free  
**Centrifuge** 1 min  
**Collect** the flow-through

**For a limited time only!**

Receive **RNase Quiet** (worth RM249.00) with purchase of any kit above.

## RNase Quiet For RNase decontamination

RNase Quiet from Nacalai Tesque is a ready-to-use solution for eliminating RNase contamination. It completely removes RNase contamination from glass and plastic equipments and laboratory tables.

## Features

- **Removes** RNase contamination effectively
- **Easy to use**, spray type
- **Easy to wipe**, no detergent required
- **Non-carcinogenic** with no DEPC



For more information, email [info@axonscientific.com](mailto:info@axonscientific.com) or call your product specialist today.



